



Product Number(s): PF30075, PF30150, PF30750

p-Fect™

Instruction Manual

Neuromics
5325 West 74th Street, Suite 8
Edina, MN 55438
Phone: 952-374-6161
Fax: 612-677-3976
Email: pshuster@neuromics1.com
Website: www.neuromics.com

Purchaser Notification

Limited License

The purchase price paid for the p-FectTM Reagent kit by end users grants them a non-transferable, non-exclusive license to use the kit and/or its separate and included components (as listed in the Kit Contents section). This kit is intended **for internal research only** by the purchaser. Such use is limited to the transfection of nucleic acid into neuronal cells as described in the product manual. Furthermore, research only use means that this kit and all of its contents are excluded, without limitation, from resale, repackaging, or use for the making or selling of any commercial product or service without the written approval of Neuromics.

Purchasers may terminate this License at any time by returning all p-Fect Transfection Reagent material and documentation to Neuromics, or by destroying all p-Fect Transfection Reagent kit components.

This document covers in full the terms of the p-Fect Reagent research only license, and does not grant any other express or implied license. The laws of Minnesota shall govern the interpretation and enforcement of the terms of this License.

Product Use Limitations

The p-Fect Transfection Reagent kit and all of its components are developed, designed, intended, and sold for research use only. They are not to be used for human diagnostic or included/used in any drug intended for human use. All care and attention should be exercised in the handling of the kit components by following appropriate research lab practices.

For more information or any comments on the terms and conditions of this License, please contact:

Neuromics
5325 West 74th Street, Suite 8
Edina, MN 55438
Email: phuster@neuromics.com

This product is manufactured for Neuromics by Genlantis, a division of Gene Therapy Systems

Contents

Introduction	3
METHODS AND PROCEDURES	3
A. DNA DILUENT SELECTION	3
B. p-Fect LIPID HYDRATION	3
C. PROTOCOL FOR USING THE DNA DILUENT	4
<input type="checkbox"/> Transfection of Adherent Cells	4
<input type="checkbox"/> Transfection of Suspension Cells	5
D. PROTOCOL FOR USING THE DNA DILUENT B	5
<input type="checkbox"/> Transfection of Suspension Cells	6
Green Fluorescent Protein	6
Secreted Alkaline Phosphatase	6
F. OPTIONAL PROTOCOL FOR LOW QUANTITY DNA TRANSFECTION	7

Kit Contents		
PF30075-75 Reactions	<i>pn-Fect Lipid Film, dried</i>	<i>1 vial</i>
	<i>Hydration Buffer, Transfection Grade</i>	<i>1 x 0.8 ml</i>
	<i>DNA Diluent</i>	<i>1 x 4 ml</i>
	<i>DNA Diluent B</i>	<i>1 x 4 ml</i>
PF30150-150 Reactions	<i>pn-Fect Lipid Film, dried</i>	<i>1 vial</i>
	<i>Hydration Buffer, Transfection Grade</i>	<i>1 x 1.6 ml</i>
	<i>DNA Diluent</i>	<i>1 x 8 ml</i>
	<i>DNA Diluent B</i>	<i>1 x 8 ml</i>
PF30750-750 Reactions	<i>pn-Fect Lipid Film, dried</i>	<i>1 vial</i>
	<i>Hydration Buffer, Transfection Grade</i>	<i>5 x 1.6 ml</i>
	<i>DNA Diluent</i>	<i>5 x 8 ml</i>
	<i>DNA Diluent B</i>	<i>5 x 8 ml</i>

p-Fect is shipped ambient and is stable stored at 4°C for up to 1 year.

Introduction

This transfection reagent is the one of most effective and widespread gene delivery reagent available. It is designed for more difficult to transfect cell types. With a choice of two optimized DNA diluent buffers, p-Fect performs effectively across a broad range of cells, offers the highest efficiencies, and is effective in the presence of serum, eliminating inconvenient media changes.

METHODS AND PROCEDURES

A. DNA DILUENT SELECTION

Select the DNA diluent suitable for your cell type as listed below; if not listed, choose the cell type that closely matches yours or consult our citations. If no data is available for your cell type, try the DNA Diluent B protocol first (Section C).

Cell Lines	DNA Diluent	DNA Diluent B	Serum
HeLa-S3	★	★★	○
HeLa	★	★★	○
COS-1	★	★	●
COS-7	★	★	●
Hep-G2	★	★	●
NIH-3T3	★	★	●
MDCK	★	★★	○
K-562	★	★★	○
CV-1	★	★	●
B15-F0	★	★	●
293	★	★	●
BHK-21	★	★	●
CHO-K1	★▲	★▲	●
PC-12	★	NR	●
P19	★	★	●
HUVEC-C	★	★	●

Legend

★ Works well	○ Works well w/o serum
★★ Works better	● Works well w/ and w/o serum
NR Not recommended	
▲ Best expression is w/o serum during 1 st hr. of transfection	

B. p-Fect LIPID HYDRATION

1. Hydrate each p-Fect lipid vial at room temperature with 0.75 ml (for T202007) or 1.5 ml (for T202015 and T202075) of the Hydration Buffer. Vortex for 10 seconds at top speed before use. Store hydrated reagent at 4°C; vortex briefly before each use.
2. Use 25 µl of the DNA Diluent or DNA Diluent B per each 1 µg of DNA. Avoid vortexing either of the DNA diluent solutions.

C. PROTOCOL FOR USING THE DNA DILUENT

3. For most cell types, use 5 μ l p-Fect per 1 μ g of DNA.
4. Your DNA can be suspended in TE buffer or purified water. A DNA concentration of at least 0.1 mg/ml works well for most reaction sizes.

- *Transfection of Adherent Cells*

5. Dilute the hydrated p-Fect reagent with serum-free medium as shown in Table 1 below.
6. Dilute the DNA with DNA Diluent as shown in Table 1 below; mix well by pipetting and incubate at room temperature for 1-5 minutes.

NOTE: Do not vortex DNA Diluent.

Table 1: Amounts of DNA, diluent, p-Fect, and Medium

DNA (μ g)	DNA Diluent (μ l)	p-Fect (μ l)	Serum-free Medium (μ l)
0.5	12.5	2.5	10.0
1.0	25.0	5.0	20.0
2.0	50.0	10.0	40.0
4.0	100.0	20.0	80.0
8.0	200.0	40.0	160.0

7. Add the diluted DNA to the diluted p-Fect reagent. Incubate at room temperature for 5 to 10 minutes to form lipid/DNA complexes (lipoplexes).

NOTE: Do not incubate for more than 30 minutes.

8. Add the p-Fect/DNA complexes directly to cells growing in serum-containing medium and incubate at 37°C. Use Table 2 below for recommended transfection volumes and DNA amounts.

NOTES: Cells plated the day before transfection should be 50%-70% confluent on transfection day. Omitting antibiotics from the media during transfection may increase expression levels; this effect is cell-type dependent and usually small. For some cells (such as HeLa S3, MDCK, CHO-K1), higher transfection efficiencies can be achieved when the initial 4-hour incubation is done in serum-free media, followed by the addition of one volume of medium containing 20% serum.

Table 2: Transfection Volumes and DNA Amounts Per Dish Size

Tissue Culture Dish Size	DNA (μ g)	Transfection Volume (ml)
96 well	0.1-0.5	0.1
24 well	0.5-2.0	0.25
6 well	2.0-6.0	1.0
60 mm	6.0-8.0	2.5
100 mm	8.0-12.0	5.0

9. Add fresh growth media as needed 24 hours post transfection; for some cell types, the old media can be replaced with fresh media at this step.

10. Depending on cell type and promoter activity, reporter gene assay can be performed 24-72 hours post transfection.

NOTE: the same protocol can be used to produce stably transfected cells: 48 to 72 hours post transfection, put the cells in fresh medium containing the appropriate selection antibiotic. Wait at least 48 hours before exposing the transfected cells to the selection medium, and for some cell types it may be necessary to wait as long as 4 to 5 days.

- **Transfection of Suspension Cells**

p-Fect reagent works well for most suspension cells, such as K562 and PC 12.

For suspension cells, the protocol is the same as described for adherent cells.

11. The day before transfection, split the cells so they are in good condition on the day of transfection.

12. While the p-Fect/DNA complexes are incubating, spin down the cells, resuspend them at 1×10^6 or 2×10^6 cells/ml in medium with or without serum, and transfer the appropriate volume to the dish as indicate in Table 3 below.

13. Add the p-Fect/DNA complexes directly to the cells, and mix well by gently pipetting 2 to 3 times.

NOTE: this step is important because some suspension cells have a tendency to clump, which reduces transfection efficiency.

14. Incubate at 37°C and proceed as described for adherent cells.

NOTE: for some hematopoietic cell lines, mitogenic agents like PHA or PMA may be added to the cells 4 hours after transfection to a final concentration of 1 µg/ml or 50 ng/ml, respectively, to enhance the levels of gene expression.

Table 3: Number of Cells and Transfection Volumes for Suspension Cells

Tissue Culture Dish Size	Number of Cells (µg)	Transfection Volume (ml)
96 well	1×10^5	0.1
24 well	0.5×10^6	0.25
6 well	2×10^6	1.0
60 mm	5×10^6	2.5
100 mm	1×10^7	5.0

D. PROTOCOL FOR USING THE DNA DILUENT B

15. For most cell types, use 3.5 µl of the p-Fect reagent per 1 µg of DNA.

- **Transfection of Adherent Cells**

16. Dilute the hydrated p-Fect reagent with serum-free medium as shown in Table 4 below.

17. Dilute the DNA with DNA Diluent B as shown in Table 4 below; mix well by pipetting and incubate at room temperature for 5 minutes.

NOTE: Do not vortex DNA Diluent B.

Table 4: Amounts of DNA, diluent, p-Fect, and Medium

DNA (μ g)	DNA Diluent B (μ l)	p-Fect (μ l)	Serum-free Medium (μ l)
0.5	12.5	2.5	10.0
1.0	25.0	5.0	20.0
2.0	50.0	10.0	40.0
4.0	100.0	20.0	80.0
8.0	200.0	40.0	160.0

18. Add the diluted DNA to the diluted p-Fect Incubate at room temperature for 5 minutes to form lipid/DNA complexes (lipoplexes).

NOTE: Do not incubate for more than 30 minutes.

19. Same as Step 8 above.

20. Same as Step 9 above.

21. Same as Step 10 above.

- *Transfection of Suspension Cells*

Same as Section ii protocol for transfection of suspension cells when using the DNA Diluent.

Green Fluorescent Protein

When green fluorescent protein (GFP) is the reporter gene used for transfections, use epifluorescence or confocal microscopy to detect expression. GFP has an excitation peak at 470 to 490 nm and emission peak at 510 nm. Expression levels of GFP can also be monitored by fluorescence-activated cell sorter analysis (FACS)³.

Secreted Alkaline Phosphatase

When heat-stable secreted alkaline phosphatase (SEAP) is the reporter gene used for transfections, use the following assay: heat supernatants from transfected cells at 65°C for 30 min to inactivate endogenous alkaline phosphatase activity. The SEAP transgene is stable during this treatment. Take aliquots of the culture media 48 hours posttransfection, and determine the SEAP activity quantitatively by using a colorimetric assay based on hydrolysis of the chromogenic substrate paranitrophenyl phosphate (PNPP). Dissolve 1 mg/ml of PNPP reagent in a solution of 1 mM MgCl₂ and 100 mM diethanolamine, pH 9.8. Add 10 μ l of 0.05% Zwittergent in PBS (free Ca²⁺ and Mg²⁺) into each well of a 96- well plate. Then add 20 μ l of the heated cell culture media to each well. For control wells, 20 μ l of water is used to normalize the volume. An alkaline phosphatase standard (EIA grade calf intestine alkaline phosphatase; Boehringer Mannheim) can be used to generate a standard curve from 10 to 10,000 pg per well. Add 200 μ l of the PNPP substrate to each well to start the enzymatic reaction. Allow the reaction to incubate at room temperature until the yellow color develops. Using 0.05% Zwittergent in PBS as the diluent virtually reduces the background to zero, which increases the detection limit of the assay. Read the plates at 405 nm using either kinetic or static mode.

F. OPTIONAL PROTOCOL FOR LOW QUANTITY DNA TRANSFECTION

The following revised protocol, 2 can be used to facilitate pipetting and transfer of DNA/lipids complexes to the cells when a low quantity of DNA ($\leq 1 \mu\text{g}$) is used for transfection.

1. Dilute hydrated p-Fect reagent with serum-free medium as in Table 5 below.
2. First dilute the DNA diluent in serum-free medium and then add the DNA. See Table 5 (A, B, and C) below for volumes of serumfree medium, DNA diluent, and DNA amounts.

Table 5: Recommended Amounts of Reagents for Optional Protocol

A: Dilution of p-Fect Reagent		
DNA (μg)	Serum-free Medium (μl)	p-Fect Reagent (μl)
0.125	49.37	0.63
0.25	48.75	1.25
0.5	47.5	2.5
1.0	45.0	5.0
B. DNA Dilution		
Serum-free Medium (μg)	DNA Diluent (μl)	DNA (μg)
46.8	3.12	0.125
43.75	6.25	0.25
37.5	12.5	0.5
25.0	25.0	1.0
C. Transfection Volume and DNA Amounts Per Dish Size		
Tissue Culture Dish Size	DNA (μg)	Transfection Volume (ml)
96 well	0.1-0.25	0.1
24 well	0.5-2.0	.25

3. Incubate 1 to 5 minutes at room temperature.
4. Proceed as in Steps 6 though 9 under the “Transfection of Adherent Cells” Section (when using the DNA Diluent).

REFERENCES

1. Felgner, JH. *et al.* (1994) Enhanced gene delivery and mechanism studies with a novel series of cationic lipid formulations. *J Biol Chem* **269**: 2550-2561.
2. Felgner, PL. *et al.* (1987) Lipofection: a highly efficient, lipid-mediated DNA-transfection procedure. *Proc Natl Acad Sci USA* **84**: 7413-7417.
3. Cheng, L. *et al.* (1996) Use of green fluorescent protein variants to monitor gene transfer and expression in mammalian cells. *Nature Biotechnology* **14**:606-609.

Do not hesitate to contact me with questions.

Pete Shuster, CEO

pshuster@neuromics.com

Direct phone: 612-801-1007